

# Bead One HF

Benchtop high frequency  
fusion instrument



**HERZOG**

# Bead One HF - Benchtop fusion instrument

## High frequency induction heating technology


The Bead One HF offers maximum flexibility and accuracy in preparation of glass beads for XRF analysis. Due to its high-frequency induction heating technology the Bead One HF enables complex multi-step fusion processes and shortens preparation time.

### Specially designed for demanding tasks


The Bead One HF offers you extensive functionality and flexibility for demanding analytical tasks. The inductive fusion method allows fast and instantaneous heating-up of the melt to the set temperature. The system operates without inertia and makes it possible to undergo complex preparation cycles requiring extremely precise temperature ramp up. This is useful for preparation procedures of, e.g., ferroalloys depending on several different oxidation and fusion steps of exactly defined temperatures and time lengths.

### Short fusion time and fast ramping up


The Bead One HF reaches temperatures of minimum 1300°C within a short period of time. Accordingly, the Bead One HF is reducing process time and increasing sample throughput. The high fusion and casting dish temperature of the Bead One HF shortens preparation time and improves bead quality for material requiring long-lasting fusion procedures. Moreover, short fusion times lead to low element losses and subsequently improve analysis accuracy and reliability. The crucible is oscillated in a circular movement to improve homogenization of the melt. Additionally, the crucible can be tilt by 45°.

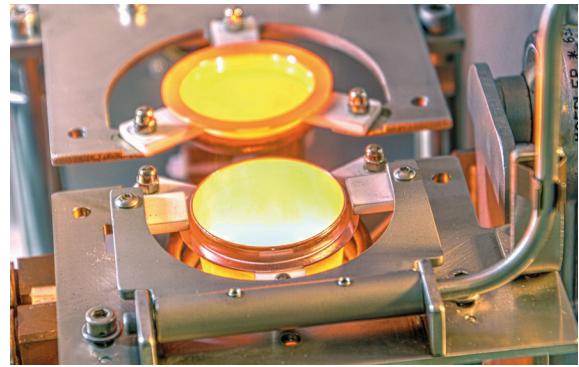

**simultaneous heating of crucible and dish**

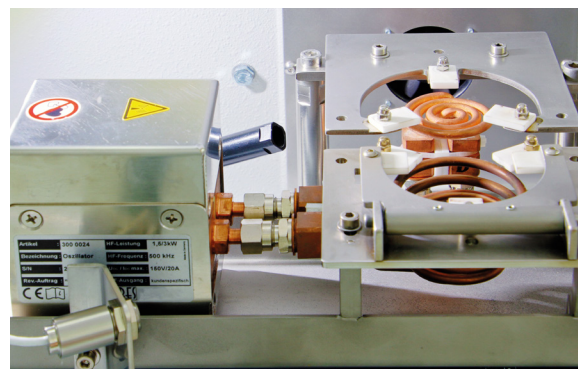


**Fusion with separate pyrometer for crucible and casting dish (optional)**

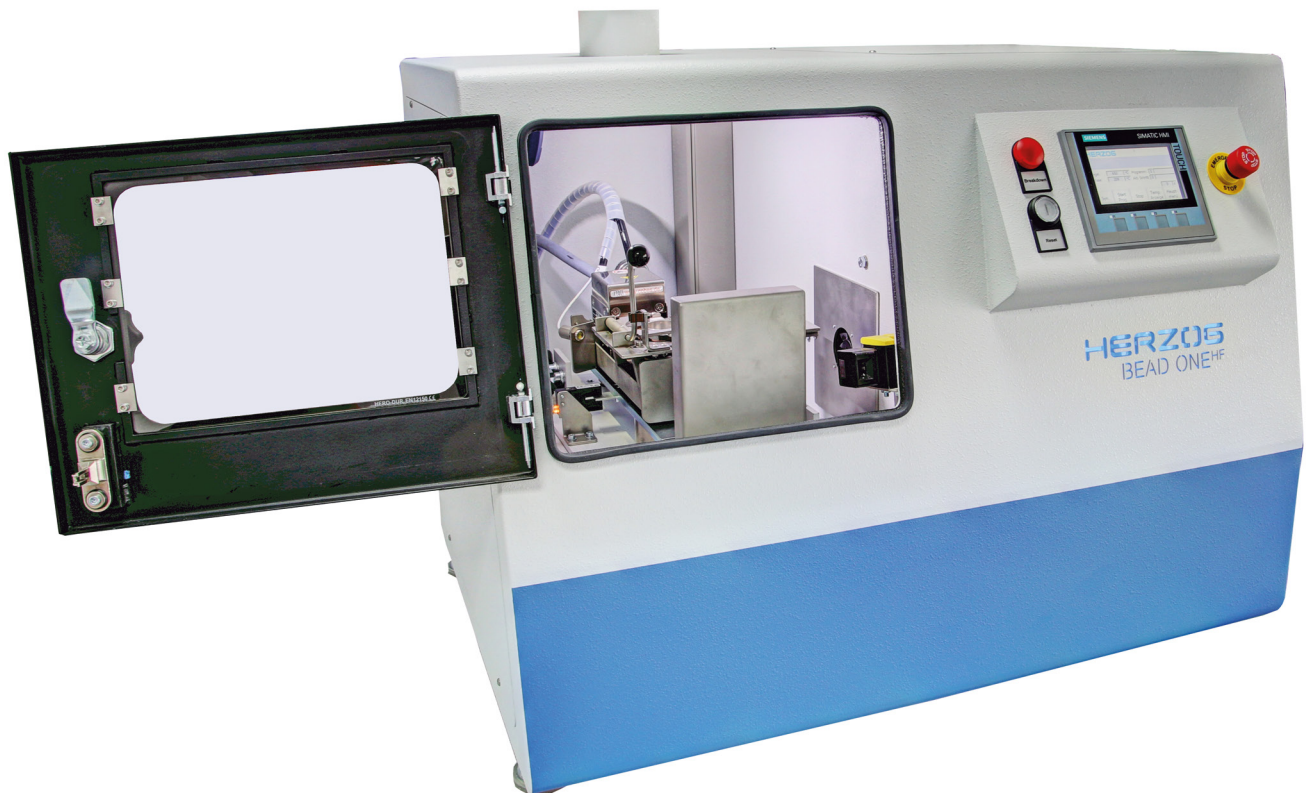

Remarkable features of the Bead One HF are the compact design, easy operation and flexible programmability. The high-frequency induction technology enables complex fusion processes within short period of time.



**Easy programm and parameter selection by HMI touch panel**

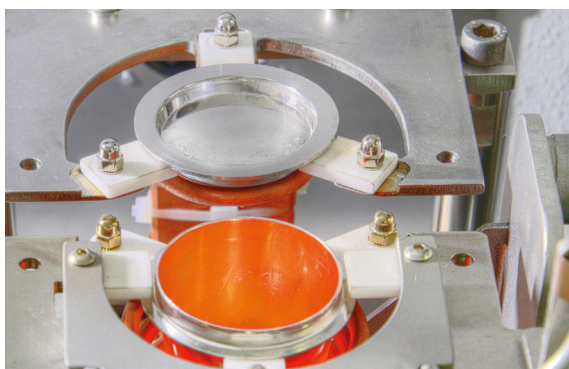

**High precision temperature control**

## Simultaneous heating of crucible and casting dish

Two separate high frequency generators supply power at the induction coil of the crucible and the casting dish. Since the casting dish is simultaneously heated up the fusion can be casted automatically into the preheated dish. Alternatively, the mold remains in the crucible until solid. The temperature curve of the dish is fully configurable and allows absolutely controlled cooling down subject to analytical requirements. Furthermore, it guarantees that glass buttons are free from reams and seeds. High performance pyrometers for crucible and casting dish ensure absolute temperature constancy. After solidifying, the casting dish is further cooled down by blowing with air.

## Safety measures

The safety door is locked during the entire fusion and cooling process. Pouring is fully automatically without need for intervention. The Bead One HF has been designed for cold-to-cold operation and the user never comes in contact with hot vessels.

## Easy to operate

Due to its small foot print design the Bead One HF can be set up flexibly at any location. Installation of the machine is very simple, no compressed air is needed. All fusion and cooling parameters can be easily set, modified, stored and monitored using the HMI panel of the machine.

## At a glance

- Benchtop fusion instrument (high frequency induction technology) for preparation of glass disks for XRF analysis
- Highly accurate and precise analysis results for a wide range of different applications
- Optimally suited for complex preparation task due to short ramping up and precise temperature control
- Easy installation and operation, full flexibility in parameter setting, and circular rocking of the crucible for perfectly fused and homogenized beads
- Maximum safety measures including safety door locking and cold-to-cold operation

# Technical data

## Bead One HF

|  |  |
|--|--|
| <b>Colour</b>                          | Blue/white<br>RAL 5007 / 7035 / 7005   |
| <b>Dimension and Weight</b>            |  |
| LxWxH                                  | 800x670x790mm  |
| Weight                                 | Approx. 140kg  |
| <b>Power supply and consumption</b>    |  |
| Voltage                                | 230 V, 50 Hz, 1/N/PE; CEE 32A  |
| Power consumption                      | 6.5 kVA  |
| Protection class                       | IP50   |
| <b>Control system and panel</b>        |  |
| Controller                             | Siemens S7   |
| Panel                                  | Touchpanel 4"  |
| Language                               | German/English   |
| <b>Fusion process</b>                  |  |
| Number of definable programmes         | 16   |
| Number of definable steps per programm | Up to 10 parameters: <ul style="list-style-type: none"><li>• Temperature</li><li>• Homogenization</li><li>• Cooling time</li></ul> |
| Temperature range crucible             | 385 – 1350°C   |
| Temperature range dish                 | 200 – 1200°C   |
| Heating of crucible and dish           | Simultaneous   |
| <b>Connections</b>                     |  |
| Cabel                                  | Terminal   |
| Compressed air                         | Not required   |
| Exhaust air stuf                       | Top of the machine; DN=80mm  |
| Cooling water; inlet/outlet            | 2x hose nozzle; NW9mm  |

## Cooling water supply

|                |   |
|----------------|---|
| Cooling water  | Min. 2l/min   |
| Inlet pressure | Min. 3.5 bar, cooling water inlet directly to the machine |
| Temperature    | 18°C – 25°C   |

## Bead diameters and cooling

|           |                         |
|-----------|-------------------------|
| Diameters | 29, 32, 34, 36 und 39mm |
| Cooling   | Electrical fan          |

## Input/output

|                      |        |
|----------------------|--------|
| Crucible, dish, bead | Manual |
|----------------------|--------|

## Options

- Temperature measuring of the dish
- Separate ultrasonic cleaning device
- Separate chiller for water



Die Ausführung der Maschine entspricht den geltenden UUV- und VDE-Vorschriften. Technische Änderungen vorbehalten.

Bead One HF/06.2017-E-1

### HERZOG Maschinenfabrik GmbH & Co. KG

Auf dem Gehren 1  
49086 Osnabrück  
Germany

☎ +49 541 9332-0  
Fax +49 541 9332-32

E-Mail [info@herzog-maschinenfabrik.de](mailto:info@herzog-maschinenfabrik.de)  
[www.herzog-maschinenfabrik.de](http://www.herzog-maschinenfabrik.de)

### HERZOG Automation Corp.

16600 Sprague Road, Suite 400  
Cleveland, Ohio 44130  
USA

☎ +1 440 891 9777  
Fax +1 440 891 9778

E-Mail [info@herzogautomation.com](mailto:info@herzogautomation.com)  
[www.herzogautomation.com](http://www.herzogautomation.com)

### HERZOG Japan Co., Ltd.

3-7, Komagome 2-chome  
Toshima-ku  
Tokio 170-0003, Japan

☎ +81 3 5907 1771  
Fax +81 3 5907 1770

E-Mail [nfo@herzog.co.jp](mailto:nfo@herzog.co.jp)  
[www.herzog.co.jp](http://www.herzog.co.jp)

### HERZOG (Shanghai) Automation Equipment Co., Ltd.

Section A2,2/F, Building 6,  
No.473, West Fute 1st Road,  
Waigaoqiao F.T.Z, Shanghai, 200131,  
P.R. China

☎ +86 21 50375915  
Fax +86 21 50375713

E-Mail [xc.zeng@herzog-automation.com.cn](mailto:xc.zeng@herzog-automation.com.cn)  
[www.herzog-automation.com.cn](http://www.herzog-automation.com.cn)

# HERZOG